LCD-Mikroskop LCD Microscope

Art. No. 15024

- DE** Bedienungsanleitung
- EN** Operating Instructions
- FR** Mode d'emploi
- NL** Handleiding
- ES** Instrucciones de uso

SERVICE AND WARRANTY:

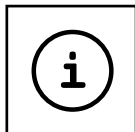
www.bresser.de/warranty_terms

MANUAL DOWNLOAD:

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
MICROSCOPE GUIDE:

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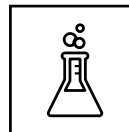EXPERIMENTS:

www.bresser.de/downloads

MICROSCOPE FAQ:

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WARNING! Not suitable for children under 3 years of age. Small parts, choking hazard. Contains functional sharp edges and points!

WARNING: Only suitable for children aged 8 years and above. Instructions for parents or other responsible persons are included and must be observed. Keep the packaging as it contains important information.

About this manual

Please read the safety instructions and user manual carefully before using the device. Keep this user manual for future reference. When selling or passing on the device, this user manual must be passed on to each subsequent owner/user of the product.

General Safety Warnings

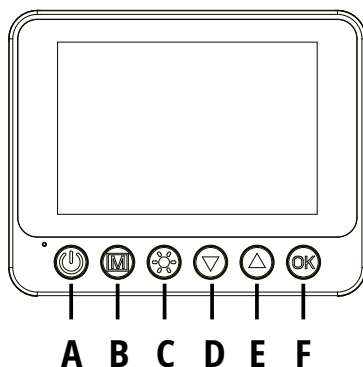
- **Choking Hazard** — Keep packaging materials such as plastic bags and rubber bands away from children! There is a risk of choking. This product contains small parts that can be swallowed by children.
- Leaking battery acid can cause chemical burns! Avoid contact of battery acid with skin, eyes, and mucous membranes. If acid comes into contact with these areas, immediately rinse with plenty of clear water and seek medical attention.
- **Risk of electric shock** — Never bend, squeeze, or pull the power and connection cables, extensions, and adapters. Protect the cables from sharp edges and heat, and inspect the device, cables, and connections for damage before use. Never use a damaged device or a device with damaged power cables. Damaged parts must be replaced immediately by an authorised service centre.
- **Risk of damage to property!** — The device must not be disassembled. In case of a defect, please contact the appropriate retailer. They will contact the service centre and may send the device for repair.
- Do not expose the device to high temperatures. Do not short-circuit or throw the device into fire! Excessive heat and improper handling can cause short circuits, fires, and even explosions!
- Do not interrupt an existing data connection to a computer until all data has been transferred. This can lead to data loss for which the manufacturer is not liable.
- The device uses a built-in 3.7V / 1000 mAh lithium-polymer battery and a 3.7V / 1500 mAh lithium-polymer battery. The batteries cannot be removed or replaced. If damaged, please contact the retailer for device repair.
- This device should only be charged under adult supervision.
- The device should be turned off while the battery is being charged.
- The product is not waterproof and should not be submerged in water or stored in a damp place.
- Press the reset button if the device freezes during operation or cannot be turned off.

Scope of Delivery




Handheld microscope, base, handle, petri dishes, box with specimens, collection box, tweezers, pipette, USB-C cable

Parts Overview


- A** On/Off switch
- B** Menu button / Rotate button / Back
- C** Adjust LED brightness
- D** Next button / Zoom out
- E** Back button / Zoom in
- F** Confirm button

- ① Screen
- ② Microphone
- ③ Charging indicator light
- ④ USB Type-C charging port 1
- ⑤ Reset button (reset device)
- ⑥ Focusing wheel
- ⑦ Adjusting the height of the microscope stage
- ⑧ Adjusting the brightness of the illuminated reticle
- ⑨ Magnetic object holder
- ⑩ Microscope stage
- ⑪ Base
- ⑫ SD card slot
- ⑬ Adjusting the height of the microscope stage
- ⑭ Adjusting the brightness of the illuminated reticle
- ⑮ Illuminated reticle
- ⑯ Speaker
- ⑰ USB Type-C charging port 2

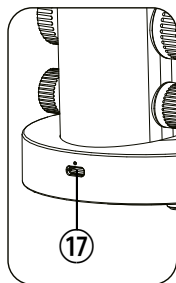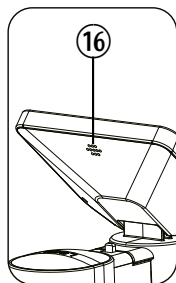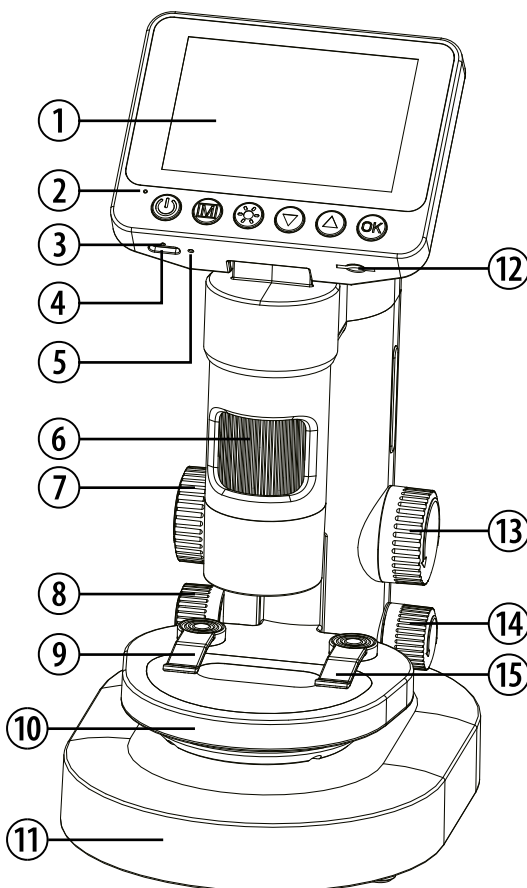Button Functions

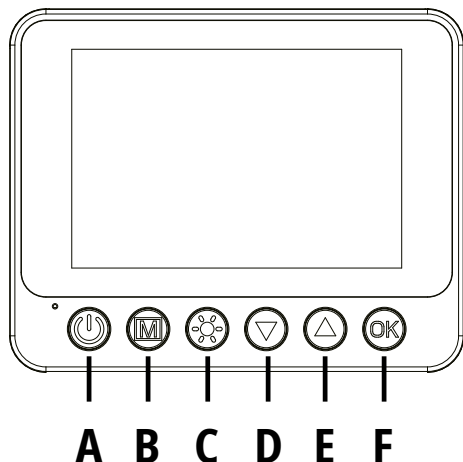(A) On/Off switch

- Press and hold the button to turn the device on or off.
- In the image gallery, briefly press the button to open the delete menu.

(B) Menu button / Rotate button / Back

- Press the menu button and use the arrow keys to choose between photo mode, video mode, playback, and settings. Press the OK button to confirm.
- Press and hold the button to vertically mirror the screen view.
- In the menu, briefly press the button to exit the menu.

(C) Adjust LED brightness

- Briefly press the button to step through the LED brightness levels (brightness 1, 2, 3, 4, 5, 6, off)

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- Briefly press the button in the menu or settings to select the next option.
- In observation mode, briefly press the button to zoom out.
- Press and hold the button to zoom faster. Zoom levels: 100X - 1000X.

(E) Back button / Zoom in

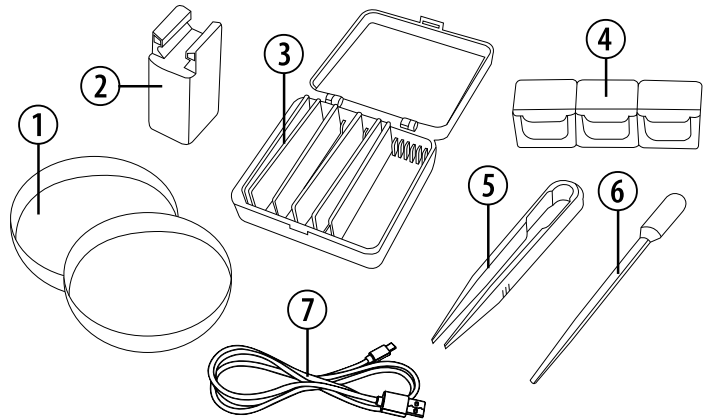
- Briefly press the button in the menu or settings to select the previous option.
- In observation mode, briefly press the button to zoom in.
- Press and hold the button to zoom faster. Zoom levels: 100X - 1000X.

(F) Confirm button



- In the menu or settings, briefly press the button to confirm the selection.
- In observation mode, briefly press the button to take a photo or record a video.

Accessories

- ① Petri dishes
- ② Handle
- ③ Box with specimens
- ④ Collection box
- ⑤ Tweezers
- ⑥ Pipette
- ⑦ USB-C cable

Handheld microscope

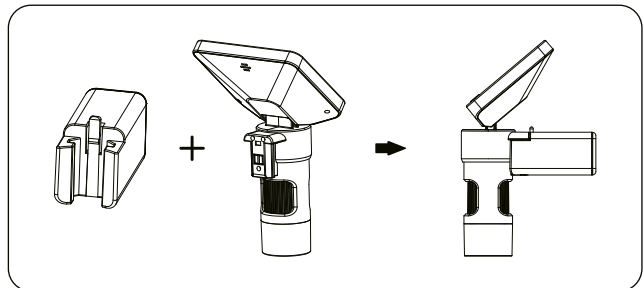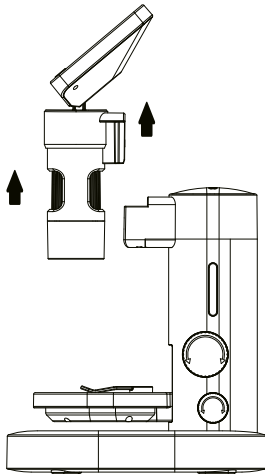
- The main part of the microscope can be removed.
- Handle and main part can be used together as a handheld microscope. Simply insert the handle into the main part.
- Tilt the screen slightly downwards and hold down the (B) menu button to vertically mirror the screen view. The microscope can now also be used as a telescope or magnifying glass.

Insert memory card

Insert the memory card (microSD/TF) in the correct orientation into the card slot until the memory card clicks into place. When removing the memory card, please press gently on the card and wait for it to pop out. You can now remove the card.

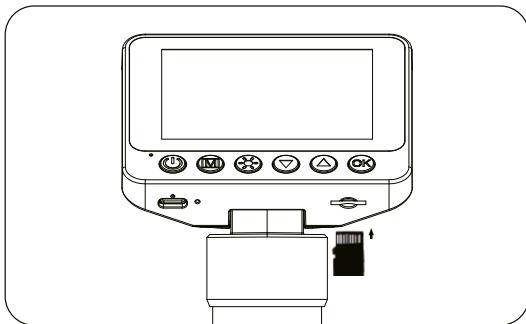
Memory cards up to 128 GB are supported.
A class 10 or higher memory card is recommended.

Do not insert or remove the card while the device is in use to avoid damage to the device or card. When you insert the memory card into the device, this card is set as the default storage device. The original information on the card may be lost. Please back up your data in advance. If the memory card is displayed as incompatible, format the card via the menu setting and try again.

Data storage and transfer

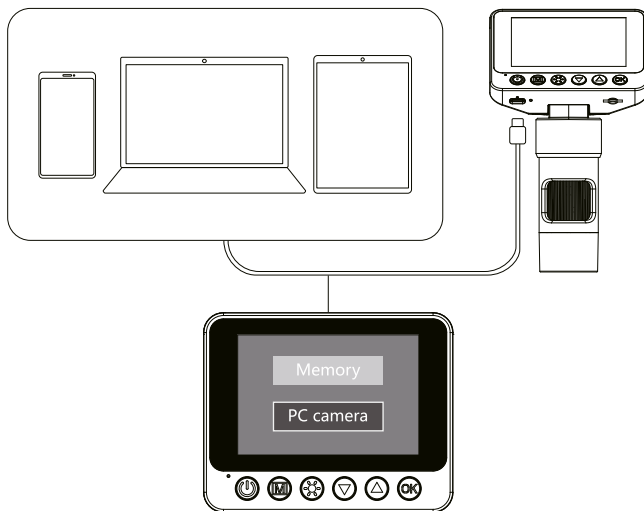
Connect the USB-C cable to a computer or smartphone to transfer data or use the microscope as a camera.

Charge the camera


Method 1:

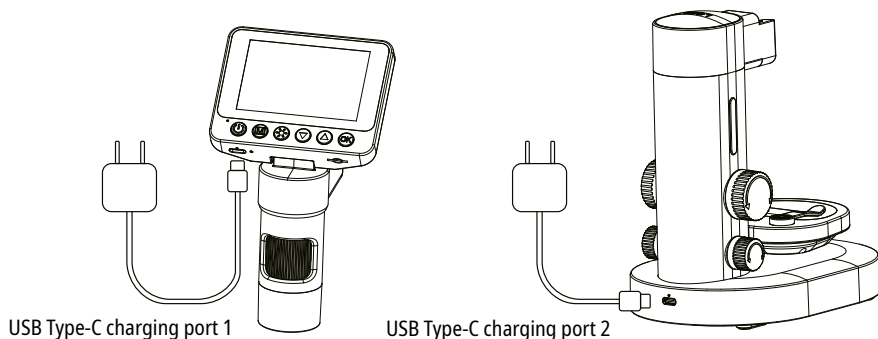Handheld microscope and base can be charged separately.

Method 2:

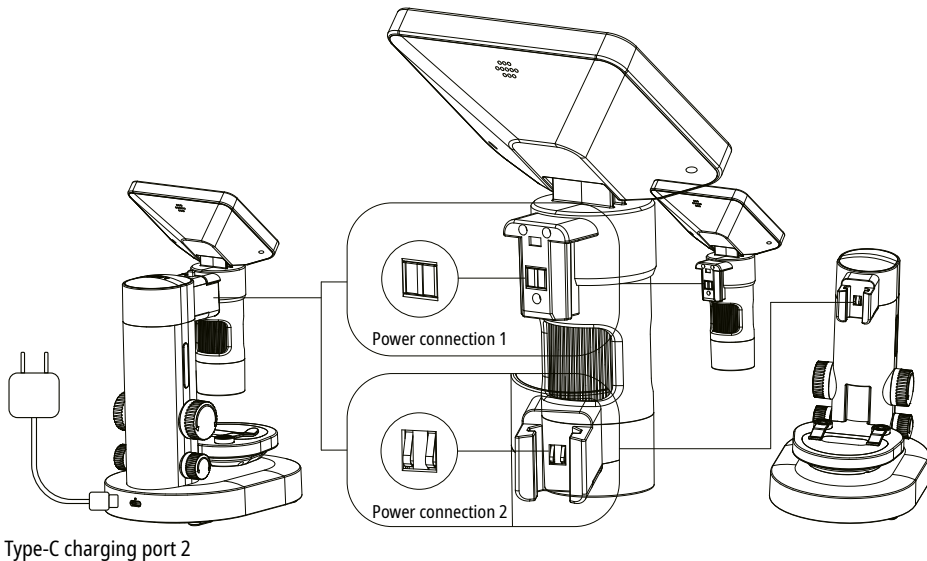When the handheld microscope and base are connected, the entire device can be charged via the USB Type-C charging port 2.

If the device's power is low, please use a 5V-1A output charger to charge the device. Using the device during charging may wear down the battery and reduce the product's lifespan. The red indicator light will be on during charging and will turn off when the battery is fully charged.

Technical Specifications	
Screen	8.9 cm (3.5") colour screen
Functions	Take photos, record videos, use as a stationary microscope, use as a handheld microscope, use as a telescope, use as a magnifying glass
Magnification	Optical Magnification: 11.4x - 64.2x Digital Magnification: 22.8x - 129.8x
Image resolution	3840x2160p / 1920x1080p
Video resolution	1920x1080p / 1280x720p
Image file format	JPG
Video file format	AVI
Dynamic frame rate	20 - 30 fps
Memory card	Max. 128 GB microSD card (not included)
Output	USB Type-C
Languages	DE, EN, CN, FR, PT, PL, NL, TR
Dimensions	120x264.5x170 mm
Weight	540 g

Experiments

You can find brochures with interesting experiments to try at the following link on the internet. How to breed prehistoric shrimp? How to make salt crystals, and much more.

EXPERIMENTS:

www.bresser.de/downloads

CLEANING TIPS

Disconnect the device from the power source before cleaning (unplug the device). Clean the exterior of the device with a dry cloth. Do not use cleaning fluids to avoid damage to electronic components. Protect the device from dust and moisture. Store the device in the supplied bag or in the original packaging.

EC Declaration of Conformity

A "Declaration of Conformity" in accordance with the applicable directives and corresponding standards has been issued by Bresser GmbH. The full text of the EC Declaration of Conformity is available at the following internet address:

www.bresser.de/download/15024/CE/15024_CE.pdf

Disposal

Dispose of packaging materials by sorting them. Please observe the current legal regulations when disposing of the device. Information on proper disposal is available from local waste disposal services or the environmental office.

Do not dispose of electrical appliances with household waste!

According to European Directive 2012/19/EU on waste electrical and electronic equipment and its implementation into national law, used electrical equipment must be collected separately and recycled in an environmentally friendly manner. Empty used batteries must be returned to battery collection points by the consumer. Further information on the disposal of devices or batteries manufactured after 6 January 2006 can be obtained from the local waste disposal service or environmental office.

Discharge the battery completely before disposal.

Warranty & Service

The regular warranty period is 2 years and begins on the day of purchase. To benefit from an extended voluntary warranty period as stated on the gift box, registration on our website is required.

The full warranty conditions and information on warranty extensions and services can be found at www.bresser.de/warranty_terms.

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Bresser UK Ltd.