Explorer Lens
Digitales Handmikroskop
Digital Hand microscope

- DE** **BEDIENUNGSANLEITUNG**
- EN** **INSTRUCTION MANUAL**
- FR** **MODE D'EMPLOI**
- NL** **HANDLEIDING**
- ES** **MANUAL DE INSTRUCCIONES**

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Hereby, Bresser GmbH declares that the equipment type with article number 15152 is in compliance with Directive: 2014/53/EU. The full text of the EU declaration of conformity

is available at the following internet address:
[http://www.bresser.de/download/15152/CE/15152_](http://www.bresser.de/download/15152/CE/15152_CE.pdf)
[CE.pdf](http://www.bresser.de/download/15152/CE/15152_CE.pdf)

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TABLE OF CONTENTS

EN

1. General Warnings	19
2. Scope of delivery	21
3. Parts overview	21
4. Operation	25
5. Preview mode (Camera mode)	26
6. Playback mode	27
7. Menu mode	27
8. Software for Windows or Mac	29
9. Cleaning and Maintenance	30
10. Disposal	30
11. EC Declaration of Conformity	31
12. Warranty & Service	31

1. GENERAL WARNINGS

- Risk of choking — Keep packaging material, like plastic bags and rubber bands, out of the reach of children, as these materials pose a choking hazard. This product contains small parts that could be swallowed by children.

- This device contains electronic components that operate via a power source. Children should only use the device under adult supervision. Only use the device as described in the manual; otherwise, you run the risk of an electric shock.
- Risk of electric shock — Never bend, pinch or pull the power and connecting cables, extensions and adapters. Protect the cables from sharp edges and heat. Before operating, check the device, cables and connections for damage. Never use a damaged unit or a unit with damaged power cables. Damaged parts must be exchanged immediately by an authorised service centre.
- Risk of property damage! — Do not disassemble the device. In the event of a defect, please contact your dealer. The dealer will contact the Service Centre and can send the device in to be repaired, if necessary.
- Do not expose the device to high temperatures. Do not short-circuit the device or throw them into a fire. Excessive heat or improper handling could trigger a short-circuit, a fire or an explosion.
- This toy contains batteries that are non-replaceable.
- Never point the light at your eyes to avoid damaging them.
- Keep instructions and packaging as they contain important information.

Check the Power supply for damage to the cord, plug, enclosure or other parts on a regular basis before use. Do not use until the damage has been repaired. The

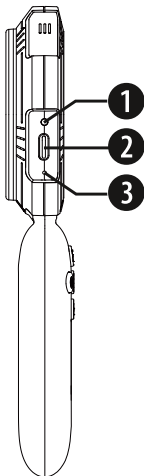
toy is not intended for children under three years. The toy must only be used with the recommended power supply. The power supply is not a toy. The toy is not to be connected to more than the recommended number of power supplies.

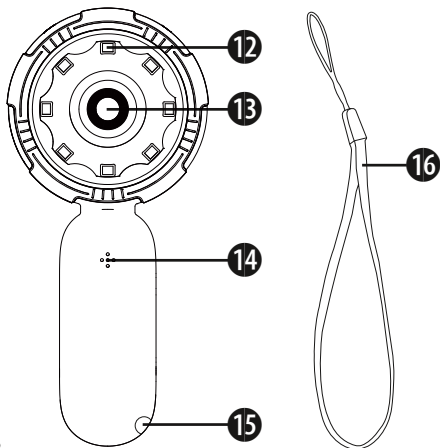
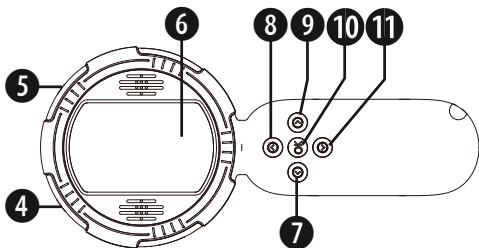
2. SCOPE OF DELIVERY

Handheld microscope, Carrying strap, USB-C cable, Instruction manual

3. PARTS OVERVIEW

1. Indicator light
2. USB Type-C connector
3. Reset button
4. Menu button
5. Power button
6. Display (5,08 cm)
7. Down button
8. Left button
9. Up button
10. OK button
11. Right button
12. LED lights
13. Camera
14. Loudspeaker
15. Hole for carrying strap
16. Carrying strap

1. **Indicator light**

When the device is powered on, the indicator light is blue. When charging, the indicator light is red. When the device is fully charged, the indicator light is green.

2. **USB Type-C port**

The USB Type-C port is provided for charging. It is also possible to connect to a PC and use the microscope as a webcam.

3. **Reset hole**

This is a small opening designed for resetting the microscope. Use a pin or a pointed object to gently press the reset button inside the hole when necessary.

4. **Menu button**

Short press to enter menu. / in Menu mode: long press to delete file

5. **Power button**

Long press for 3 seconds to turn the device on or off.

6. **Display 5,08 cm (2")**

This is the central area for you to view, configure and execute the microscope's capturing functions. The images transmitted by the objective lens will be displayed on the screen for your observation. You can magnify the images by zooming/ out for detailed viewing and capturing.

7. **Down button**

In preview mode: Zoom out / in Menu mode: down button

8. **Left button**

In preview mode: LED brightness down / in Menu mode: left button

9. **Up button**

In preview mode: Zoom in / in Menu mode: up button

10. **OK button**

In preview mode: short press to take a picture / in Menu mode: confirmation button

11. **Right button**

In preview mode: LED brightness up / in Menu mode: right button

12. **LED lights**

If there is not enough light in the observation environment, you can increase the brightness. Use the right or left button to increase or decrease the brightness.

13. **Camera**

14. **Loudspeaker**

4. OPERATION

Charging:

Insert the USB Type-C plug into the Type-C port of the microscope, and connect the other end to a USB power supply for charging (Power supply not included) or a Computer. The charging time is about 2 hours, and the usage time is about 5 hours. The charging indicator will be red while charging and will be green when fully charged.

Port: USB-Type-C (5V/1A)

Power consumption: 200mA, 1W

Battery: 650mAh (3.7V)

Battery manufacturer:

Shenzhen Yong Jia Neng Technology Co., Ltd, East of 3rd Floor, Building 3, Wangye Industrial Plant, Baotian 3rd Road, Xixiang, Baoan District, Shenzhen, Guangdong, China. Mail: 542258801@qq.com
.....

Power on:

Press and hold the On/off button for 3 seconds. The start screen appears and after a while you will be taken to the preview mode.

*Note: If you connect the USB-C cable to a PC you will have the option to use the microscope as "PC Camera" or to enter "Charging Mode". **PC Camera** will enable you to use the microscope with the "HiView" Software available for Windows or Mac. **Charging mode** allows you to use the camera while it is charging.*

Menu navigation:

Press the menu button to enter the menu mode. Press left or right button to navigate through the menu. Press the OK button to confirm your selection.

5. PREVIEW MODE (CAMERA MODE)

LED Illumination:

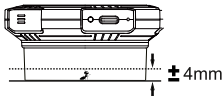
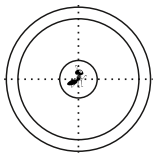
When in preview mode, Press the right or left button, to turn the brightness up or down.

Magnification:

When in preview mode, press the up or down button to zoom in or out.

Taking photos:

When in preview mode, aim the microscope at the object to be observed, make sure the object is centered under the translucent cover. The device has an integrated 4 MB flash memory for storing your images.

6. PLAYBACK MODE

Photo Playback:

When in preview mode, press the Menu button to enter the menu. Use the left button to navigate to "Playback". Press the OK button to confirm. You are now in the playback mode.

Delete files:

While viewing a file in playback mode, long press the menu button to access the file delete option.

Return to the menu:

While in playback mode, short press the menu button to return to the menu.

7. MENU MODE

Playback:

View your captured images, which are stored on the 4 MB flash memory.

Language:

Set the language: German, English, French, Dutch, Spanish

Brightness:

8 levels of brightness can be selected.

Auto Power Off:

After selecting the automatic shutdown time, the microscope is automatically switched off when it is not in use. If automatic shutdown is not required, please select 'Off'.

Screensaver:

Select the time until the LCD screen switches off automatically to save battery life.

Volume:

Adjust the volume of the loudspeaker. Select "0" to mute.

Format:

Delete all user files saved in memory.

Default Settings:

Reset to factory default settings.

Version:

Check firmware version.

8. SOFTWARE FOR WINDOWS OR MAC

EN

Software download:

Please download the applicable software for your system.

HIVIEW SOFTWARE DOWNLOAD:

www.bresser.de/download/15152
.....

You can also search for “HiView” in the Microsoft Store or App Store.

Alternative Link: <http://soft.hvscam.com>
.....

How to use:

- Step 1: Connect the microscope to the computer via a Type-C cable.
- Step 2: Start the HiView Software.
- Step 3: Long press the microscope power button to turn the device on.
- Step 4: The microscope gives you the option to choose between “PC Camera” and “Charging mode”. Choose “PC Camera” to use the microscope as a webcam.
- Step 5: In the HiView Software under “Device” choose “Digital Microscope”
- Step 6: Click on “Photo” to take a photo or “Video” to capture a video.

Note: While in this mode, only the right and left buttons are functional (to turn the brightness up or down). Long press the power button to shutdown the device.

9. CLEANING AND MAINTENANCE

- Before cleaning the device, disconnect it from the power supply!
- Only use a dry cloth to clean the exterior of the device. To avoid damaging the electronics, do not use any cleaning fluid.

10. DISPOSAL

Dispose of the packaging materials properly, according to their type, such as paper or cardboard. Contact your local waste-disposal service or environmental authority for information on the proper disposal.

Do not dispose of electronic devices in the household garbage!

As per Directive 2002/96/EC of the European Parliament on waste electrical and electronic equipment and its adaptation into German law, used electronic devices must be collected separately and recycled in an environmentally friendly manner. Empty, old batteries must be disposed of at battery collection points by the consumer. You can find out more information about the disposal of devices

or batteries produced after 6 January 2006 from your local waste-disposal service or environmental authority.

Discharge the battery completely before disposal.

11. EC DECLARATION OF CONFORMITY

Hereby, Bresser GmbH declares that the equipment type with article number 15152 is in compliance with Directive: 2014/53/EU. The full text of the EU declaration of conformity is available at the following internet address:
http://www.bresser.de/download/15152/CE/15152_CE.pdf

12. WARRANTY & SERVICE

The regular warranty period is 2 years and begins on the day of purchase. To benefit from an extended voluntary warranty period as stated on the gift box, registration on our website is required.

You can consult the full warranty terms as well as information on extending the warranty period and details of our services at

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BRESSER Benelux B.V.
Bresser UK Ltd.