BRESSER® **MM-23**

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
GB INSTRUCTION MANUAL
NL HANDLEIDING
DE BEDIENUNGSANLEITUNG

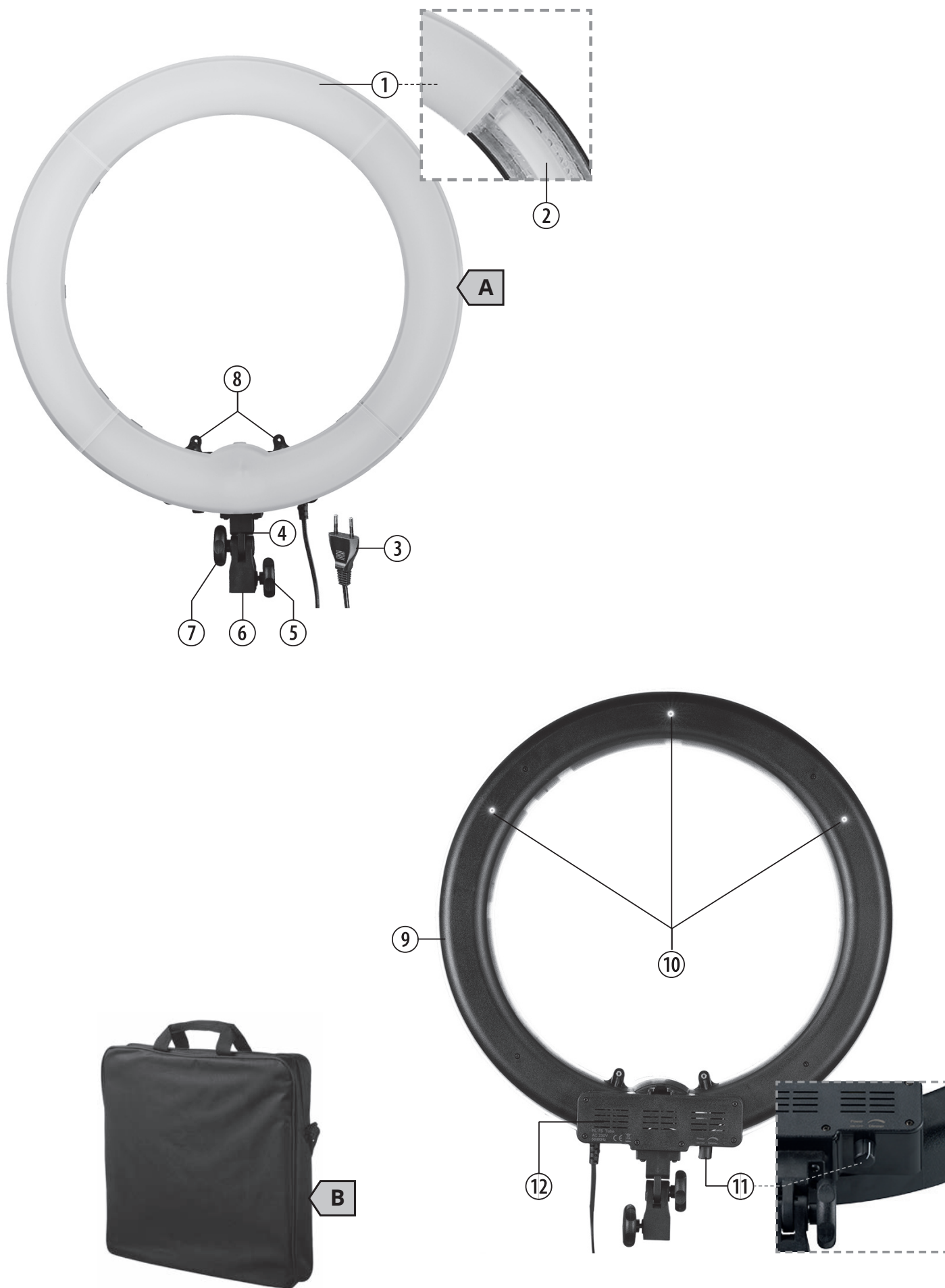Fig. 1

**GB INSTRUCTION MANUAL****GENERAL WARNINGS****⚠ DANGER OF BODILY INJURY!**

Do not use in the presence of children or animals. Store out of reach of children and animals.

Allow the unit to cool down for some time before touching it. During operation, individual components may generate a lot of heat. Touching it can then lead to painful skin burns!

⚠ RISK OF MATERIAL DAMAGE!

Make sure that no one can trip over connection cables of the device. The device could be knocked over.

Do not expose the device to moisture. Moisture penetrating the interior can cause irreparable damage to the electronics.

Do not subject the device to excessive vibrations.

The manufacturer accepts no liability for damage to property caused by improper or insufficiently safe installation!

⚠ DANGER OF AN ELECTRIC SHOCK!

This device contains electronic parts that are powered by a power source (power adapter). Only use the device as described in the manual. Otherwise, you run the DANGER of an ELECTRIC SHOCK.

Never bend, squeeze or pull power cables, connecting cables, extensions or connectors. Protect cables from sharp edges and heat. Check this device, cables and connections for damage before use. Never attempt to operate a damaged device, or a device with damaged electrical parts! Damaged parts must be replaced immediately by an authorized service agent.

Operate the device only in a completely dry environment and do not touch the device with wet or damp body parts.

⚠ RISK OF GLARE!

Make sure that people and animals do not look into the light. The high luminosity can cause glare to the eyes and afterimages. This impairment of vision can lead to accidents.

⚠ FIRE/EXPLOSION HAZARD!

Always place the device in such a way that it cannot tip over and ensure that it is stored in a reasonably safe place.

• Keep children and animals away from the device! They could knock the device over.

- Make sure that no one can trip over connection cables. The device could be knocked over.
- Do not expose the device to high temperatures. Do not short-circuit the device or dispose of in fire! Excessive heat and improper handling can cause short circuits, fires and even explosions!
- Never cover ventilation slots - not even for a short time! Heat build-up can cause fires.

! IMPORTANT NOTE!

Do not disassemble the device! In the event of a defect, please contact your dealer. They will contact the Service Center and can arrange the return of this device for repair if necessary.

Disconnect the power supply by pulling the mains plug when the appliance is not in use, in case of a longer interruption of operation and before any maintenance and cleaning work.

Place your device so that it can be disconnected from the power supply at any time. The power outlet should always be near your appliance and should be easily accessible, as the plug of the power cord serves as a disconnect device from the mains supply. To disconnect the unit from the mains, always pull the mains plug and never pull the cable!

**WARRANTY**

www.bresser.de/warranty_terms

PARTS OVERVIEW (Fig. 1)

- ❶ Semi-transparent cover (four-part, removable) ❷ Illuminant, ring-shaped¹
❸ Euro power plug ❹ Tilt head ❺ Knurled screw (tripod connection) ❻
Tripod connection ❼ Knurled screw (tilt head)
❽ Connection preparation for camera mount* ❾ Lamp housing ❿
Function lights ⓫ On/Off/Dimmerknob ⓬ Ventilation slots

SCOPE OF DELIVERY (Fig. 1)

Ring light with tilt head and power cord/plug (1x)(A), bag (1x)(B)

PUTTING THE RING LIGHT INTO OPERATION AND USING THE DIMMER

1. Insert the plug of the mains power cable into the wall socket.
2. Turn the on/off/dimmer knob (11) to the right until a slight click is heard. The light is on and glowing dimly.
3. Turn the on/off/dimmer knob to the right or left to continuously adjust the desired light intensity.

NOTE: Turning the control knob to the right increases the light intensity, turning it to the left decreases the light intensity. Depending on the light intensity set, one, two or three function lights come on.

4. To turn off the ring light, turn the on/off/dimmer control knob to the left until you hear a slight click. The lamp goes out.

MOUNTING THE RING LIGHT ON A TRIPOD*

IMPORTANT: Hold the ring light firmly during installation to prevent damage from falling.

NOTE: A suitable tripod* with spigot attachment* is required to use the ring light on a tripod.

1. Unscrew the knurled screw on the tripod connection (5) as far as possible so that the tripod connection (6) is exposed.
2. Place the ring light with the sativa connection (6) on the spigot attachment* of the tripod*.
3. Hand-tighten the knurled screw (5) again to securely mount the light on the stand* without wobbling.

TILTING THE RING LIGHT

IMPORTANT: Hold the ring light firmly while tilting to avoid damage from sudden flipping.

1. Loosen the knurled screw on the tilt head (7) so that the ring light can be tilted over the tilt head (4).
2. Move the ring light to the desired position and hand-tighten the knurled screw (7) again to securely fix the light.

CLEANING

Disconnect the power supply before cleaning! (switch off the device and pull out the mains plug)

Clean the device only on the outside with a dry cloth. To avoid damage to the electronics do not use cleaning fluids.

EC DECLARATION OF CONFORMITY

CE A "Declaration of conformity" in accordance with the applicable directives and corresponding standards has been prepared by Bresser GmbH. This can be viewed at any time upon request at www.bresser.de/download/F001428/CE/F001428_CE.pdf

UKCA DECLARATION OF CONFORMITY

CE A "declaration of conformity" in accordance with the applicable directives and corresponding standards has been prepared by Bresser UK Ltd. This can be viewed at any time at www.bresser.de/download/F001428/UKCA/F001428_UKCA.pdf

Bresser UK Ltd., Suite 3g, Eden House, The Industrial Estate, Edenbridge TN8 6HF, United Kingdom

TECHNICAL DATA

Model	F001428
Type	fluorescence ring light, dimmable
Housing colour	black
LED illuminant	fluorescence light tube, ring-shaped ¹
Power	75W ²
Light colour	5,500 K
Light output	approx. 300 lm
Power supply	220V, via mains cable/plug
Material	metal/plastic
Dimensions	inside diam: 37.5 cm outside diam.: 48.0 cm

¹The light source is permanently installed at the factory.

²Comparable to 375W for a conventional halogen lamp.

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GB IE

Please contact the service centre first for any questions regarding the product or claims, preferably by e-mail.

e-mail: service@bresseruk.com
Telephone*: +44 1342 837 098

BRESSER UK Ltd
Customer Support
Suite G3, Eden House
Enterprise Way
Edenbridge, Kent TN8 6HF
United Kingdom

*Number charged at local rates in the UK (the amount you will be charged per phone call will depend on the tariff of your phone provider); calls from abroad will involve higher costs.

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BRESSER Benelux B.V.